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- *Biocide with sporicidal activity*
- *Activity proved in short time*
- *Simplicity of use*
- *Low odour*

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INDICATIONS

Disinfection of floors, walls and equipment in the Pharmaceutical Industries for clean rooms (A and B grades).

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Biocidal activity on field surfaces and strains according to EN 13697.

Other data concerning the product (material compatibilities...): see scientific file.

For each batch, Certificate of gamma irradiation of packaging items, Certificate of analysis, LAL Testing analysis and Sterility report are available.

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- pH: 2.3
- Density: 1.0473
- Refraction index: 1.3456

INSTRUCTIONS FOR USE

Concentrated product to dilute in the ANIOS DW IP STERILE* dilution water can.

Used on previously cleaned surfaces, at a dilution of 5%. Remove the first package before entering into the clean room. Remove the second package on first use. Proceed with the dilution.

For 5L solution :

For the totality of the **250ml bottle in the ANIOS DW IP STERILE* 4.75L can.**

For 1L solution :

For the totality of the **45ml bottle in the ANIOS DW IP STERILE* 855ml bottle.**

Homogenize the solution.

Evenly apply over the surface to be treated with a suitable cleaning wipe, in sufficient amount (+/- 30 ml/m²). Ensure that the surfaces remain wet throughout. Respect the indicated contact time for the required antimicrobial activity. If necessary remove the surplus after the contact time. For frequency of use and cleaning of the application equipment, please refer to hygiene plan in place. Any preparation not used within 8 hours (open bottle) or 5 days (closed bottle) after dilution will be disposed of. Write the information regarding dilution on the ANIOS DW IP STERILE⁽¹⁾ label. At the dilution of use, the product doesn't show incompatibility with commonly found materials.

PRECAUTIONS FOR USE

Dangerous - respect the precautions for use (drawn up according to the European rules in force regarding the classification and labelling of chemical products). Use the biocides with precaution. Before use, read the label and the product information.

Store upright between +5°C and +25°C.

Biocidal product used for surface disinfection (Group 1 - PT 2) - reserved for professional use.

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